

# Androiscope

AM150X / AM200X / AM400X



[\[ App. Download \] – down.unimakes.com](http://down.unimakes.com)

or

Open “Play Store”, Search “DroidView”, Download & Install it

## Contents

---

1. Microscope x1
2. Small metal Stand x1
3. Micro-Ruler x1
4. Product Guide x1
5. OTG Adapter x1

## Caution

---

-  This product supports the Android devices of version 3.1 or above, and the devices must support OTG function.
-  The OTG function must be tested with a U-disk, a mouse and a keyboard via an OTG adapter, they must work very well before using to this Androiscope. Otherwise, the OTG function is not qualified for the Androiscope. Please change another Android device.
-  Please use the enclosed OTG Adapter and plug correctly until the end of the OTG Adapter.
  -  Connections can cause damage to the device, the screen is turned off or broken images.
-  Connect the cable is not pulled tight.
  -  Connections are disconnected or cable may be damaged.
-  Do not drop the product on the floor or use excessive shock.
  -  It can cause damage to the lens.
-  After use, please put the protective cap.
  -  Dust can make difficult to observe objects.

## How to use

---

### STEP 1. Download and Install Application

- Scan the QR CODE and download the "DroidView" APP and install it; you can also download it from the Google Play Store or directly visit our website [down.unimakes.com](http://down.unimakes.com) to download it.



### STEP 2. Connect to Android Devices

- Attach the enclosed OTG gender.
- Connect the Androiscope on Android-based mobile devices (Smartphone, Tablet PC, etc.).
- Application will be executed automatically.

### STEP 3. Image Preview & Capture

- Approach the microscope close to the observation target.
- Adjust the focus of the object using the wheel.
- Fit the image by moving the focus a little bit if the image is not clear.
- Press capture button if you want to capture the image.

### Product Warranty

---

We conduct quality assurance and follow-up services in accordance with regulations prescribed in the consumer damages and Consumer Protection..

If, when a manufacturing defect or a fault arises spontaneously within the warranty period will be exchanged through the dealer for repairs or parts or customer support center A/S center.

1. This product is made through a strict quality control and inspection.
2. You can get the organic free/non-free A/S in the event of one year after the purchase of this product. However, in the case of capital A/S, even within the warranty period, you will need to pay a fee according to the service history (service fee includes repairs, component costs, travel expenses and other additional costs.).
3. This warranty is only for domestic use and, not be re-issued to well-kept with manual.
4. If you require a repair, you should by all means bring and submit a certificate.
5. If a malfunction occurs in the product, please contact the Customer Support A/S center.

### Non-Free service case

---

1. The product warranty period ends
2. Consumer intentional, in the event of a failure due to negligence (e.g., foreign substance inputs, directly or indirectly external shocks)
3. In the event of any modifications due to failure (for example, the decomposition of the product, modifications, customization options, and unused consumables)
4. In the event of a failure due to unstable power supply (e.g., power supply phase, connected equipment failure)
5. In the event of failure due to natural disasters (e.g., fire, salt, gas, earthquake, storm and flood damage)

☞ The service material retention period is three years after we stop selling the product.